

How to Compare Two Images On a Mac

1. Connect the SmartMicroScope to the USB port and then open the software.


2. Using the Tool Menu on the right Click on the Compare Tool



3. Two windows will open on the desktop

4. Select the images that you would like to compare from the "File Browser"

5. The First image will appear on the left in the SmartMicroScope Viewer for Mac

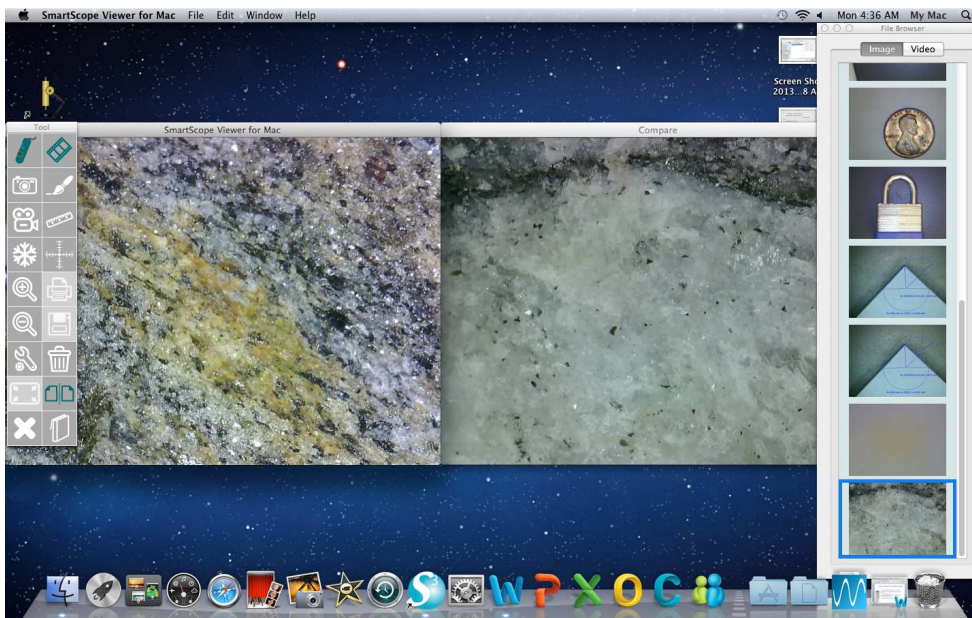

6. Click on the second image and double click on the Compare window to insert the image

7. Click the Save Tool



to save to file or Click on the Print Tool to print



8. Click the Compare Tool again to return to one screen.

*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com

1-866-722-2051

 SMARTSCHOOL
s y s t e m s

www.smartschoolsystems.com