


How to Measure A Triangle Using the SmartMicroScope 5M on Mac

Connect the SmartMicroScope 5M to the USB port and then open the software


Place the SmartMicroScope 5M on a ruler and adjust the focus knob until the image is sharp.



1. Using the function keys at the side of the Viewer, Click on the Measurement Tool
2. From here, The Measurement Window will Open

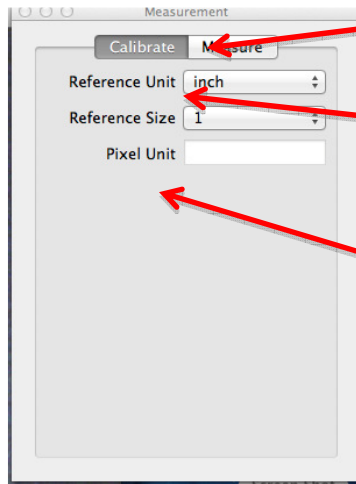


3. Click the Calibrate Tool
4. Change the Reference Unit to appropriate measurement (inch, mm, mil)
5. Change the Reference Size to 10, depending on the measurement

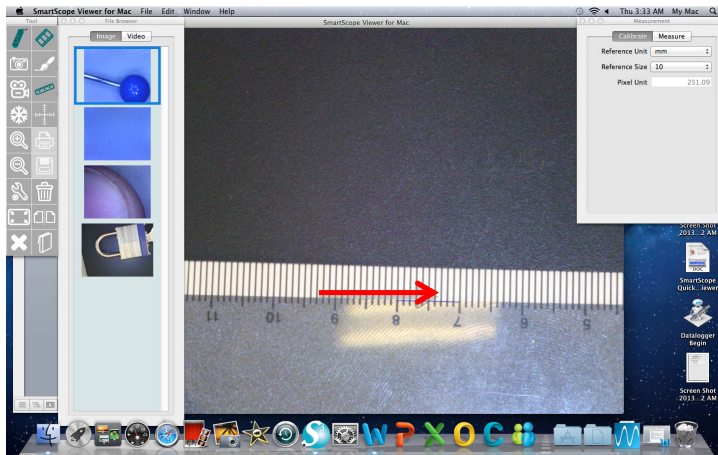


6. Using the mouse click on the Left side of the ruler and drag to the right
7. Release the mouse and your calibration is complete

*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com


How to Measure A Triangle Using the SmartMicroScope 5M on Mac

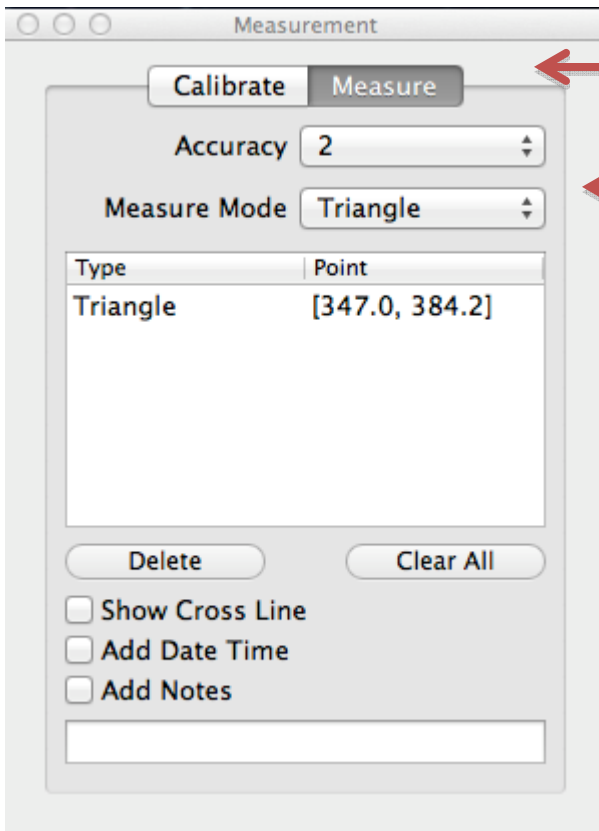


8. Click on the Measure Tool
9. Choose the Measure Mode

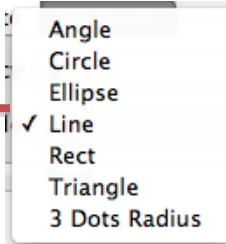

- 10.
11. Place your image under the SmartMicroScope

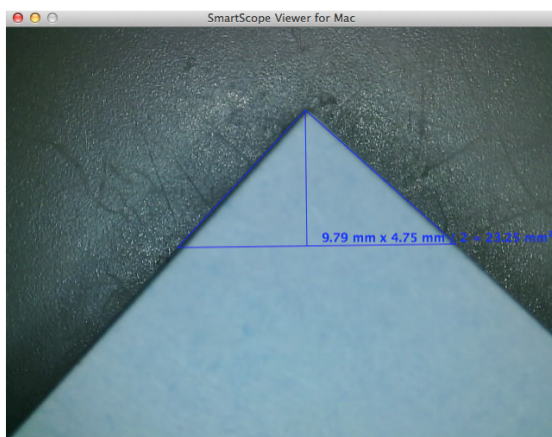

12. Place your mouse at the top of the angle
13. Click and hold the mouse and drag down the side of the image
14. Release the mouse and drag the mouse across the image
15. Click and Release the mouse and the measurement will be displayed

16. After finishing the measurement click the measurement Tool  to return to viewer

*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com