


How to Measure a Line Using the SmartMicroScope 5M on Mac

Connect the SmartMicroScope 5M to the USB port and then open the software


Place the SmartMicroScope 5M on a ruler and adjust the focus knob until the image is sharp.



1. Using the function keys at the side of the Viewer, Click on the Measurement Tool
2. From here, The Measurement Window will Open

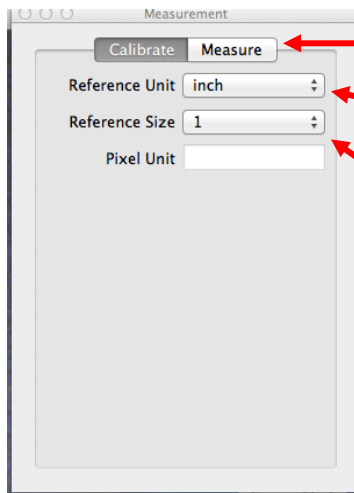


3. Click the Calibrate Tool
4. Change the Reference Unit to appropriate measurement (inch, mm, mil)
5. Change the Reference Size to 10, depending on the measurement
6. Place your ruler under the SmartMicroScope bring into focus

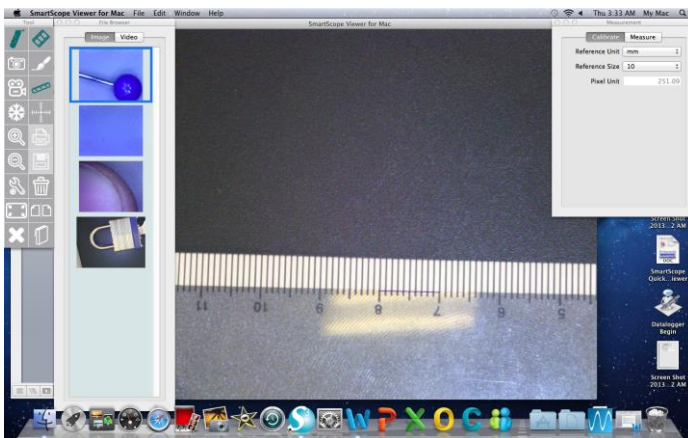


7. Place your ruler under the SmartMicroScope
8. Using the Focus Dial bring into focus
9. Using the mouse click on the Left side of the ruler and drag to the right
10. Release the mouse and your calibration is complete

****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com

How to Measure a Line Using the SmartMicroScope 5M on Mac

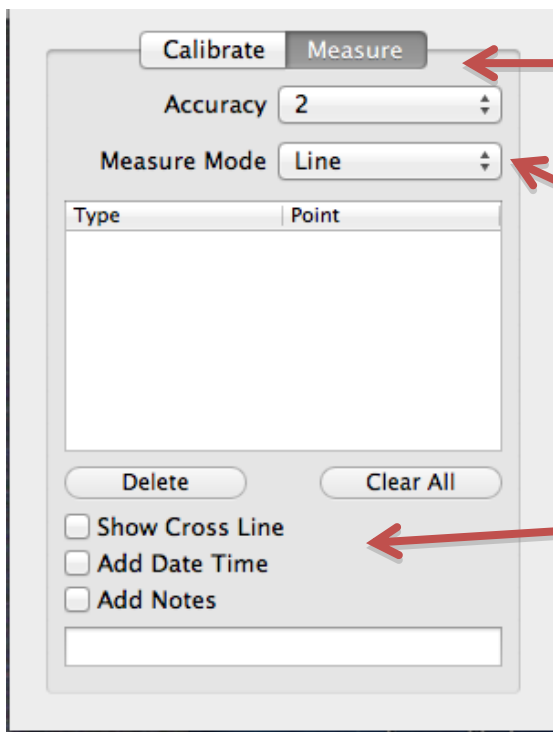


11. Click on the Measure Tool
12. Choose the Measure Mode

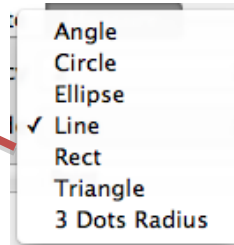

12. Click on the "Show Cross Line (when needed)

13. Now begin measuring your image

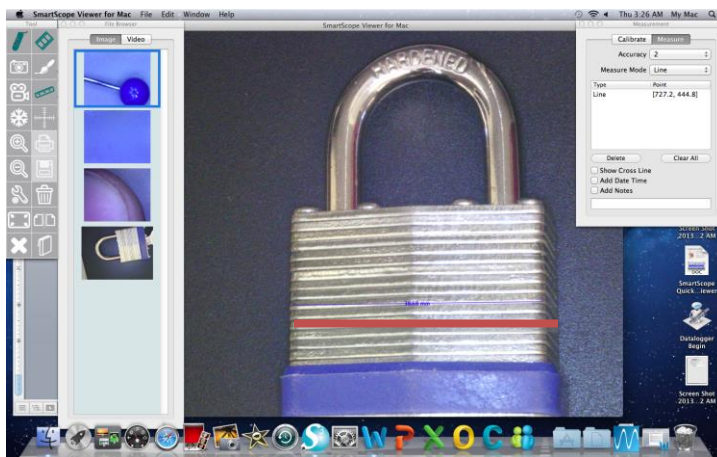

14. Move the cursor to where you want to begin measuring

15. Click and hold the mouse


16. Drag the line across the image

17. Release the mouse and the measurement will be displayed

To Save your image click the Save Tool



or Print with the Print tool

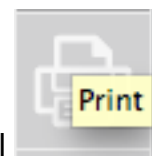

18. After finishing the measurement click the Measurement Tool



to return to viewer

*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com