

How to Compare Two Images on the SmartMicroScope 5M

1. Connect the SmartMicroScope to the USB port and then open the software.
2. Using the function keys at the top of the SmartMicroScope Viewer 5M Click on the Compare Tool



3. Two windows will open in the viewer
4. Select the images that you would like to compare from the bottom of the screen



5. The first image will appear on the left
6. Drag and drop the second image in the right window
7. Click the Save button to save for later or



8. Click on the Compare Tool to return to a single image screen

****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com

How to Compare Two Images on the SmartMicroScope 5M

Compare Two Live Images


(Requires two SmartScope 5Ms)

1. Connect both SmartMicroScopes to the USB port and then open the software. (The software must be opened with both scopes plugged in)



2. Click on the Compare Tool

3. Two screens will appear

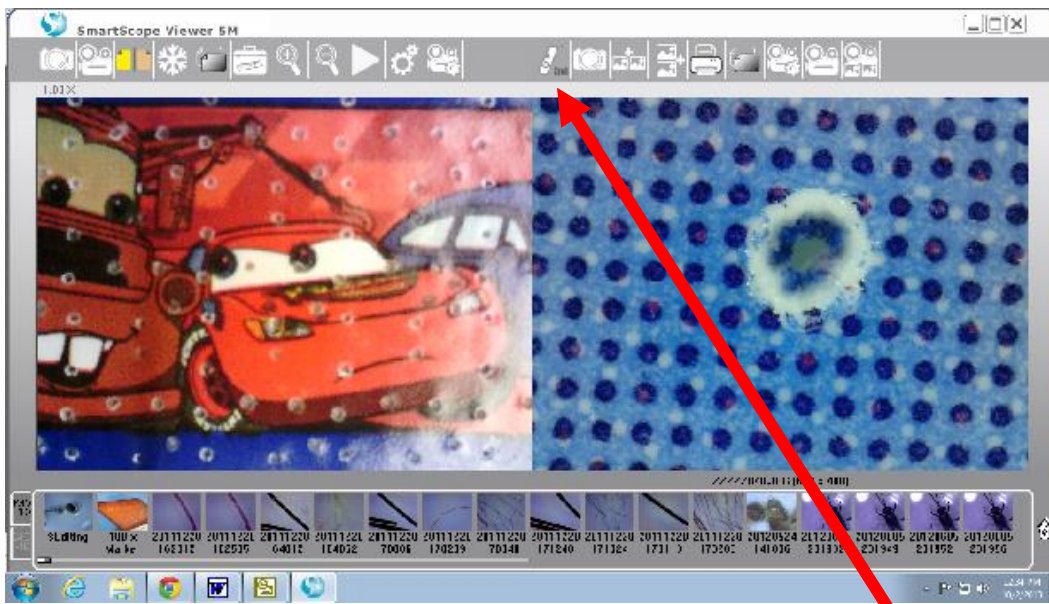

4. A Second tool bar will appear, Click on the 2nd Image Tool
5. Turn the focus dial from each scope to bring them both into focus. From here you can take a picture or record videos from both scopes at the same time.
6. When finished press the Compare Tool again to return to one screen

****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com