


# How to Create Time Lapse Video Using the SmartMicroScope 5M

1. Connect the SmartMicroScope to the USB port and then open the software.
2. Using the function keys at the top of the SmartMicroScope Viewer 5M Click on the




1. The Video Setup Window will open

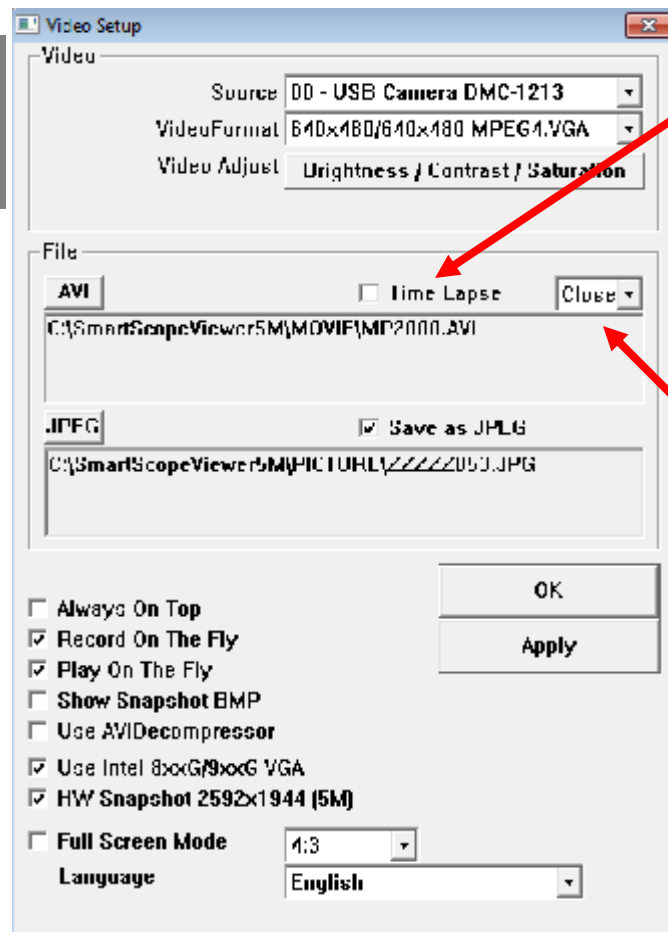

2. Click the Time Lapse tool

3. Click on the Close drop down box and choose your setting

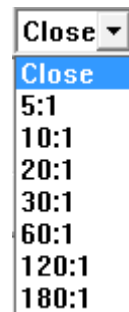

4. Press OK and your changes will automatically be added to the screen
5. Beginning recording your time lapse by pressing the Video Record icon on the toolbar

**NOTE:** If you choose 5:1, you must let the video record for at least 5 seconds. If 180:1 is chosen, it must run for at least 180 seconds to store the video.

\*\*\*\*\*Technical support, please call 281-312-1297, 1-866-722-2051 or email [support@smartschoolsystems.com](mailto:support@smartschoolsystems.com)