


Where Are My Images and Videos?

Connect the SmartMicroScope to the USB port and then open the software.


1. Using the function keys at the top of the SmartMicroScope Viewer 5M, Click on the Video Setup Tool



2. From here, The Video Setup window will open

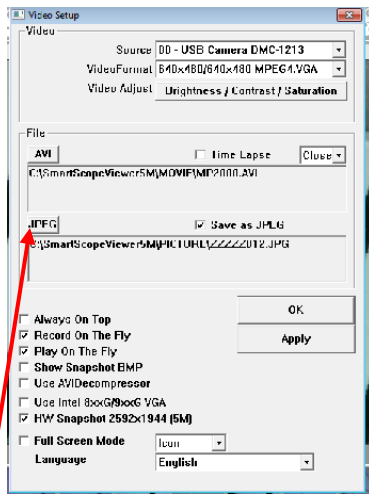

3. You will see where both your videos and images are stored.
4. If you want to change where the images are stored, Click on the JPEG icon and the “Save As Window” will open.

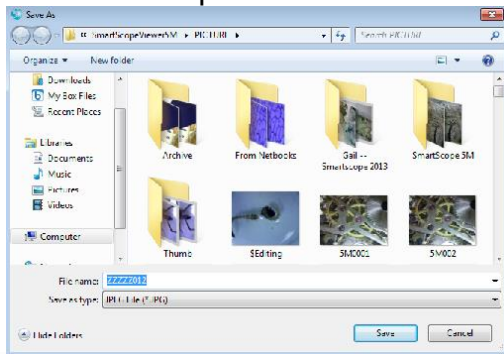

5. You can now redirect the software to store your images in a new location and change how the images are named.

NOTE: Videos can be located and redirected in this same way.

*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com