

# SmartMicroScope iGo2 Quick Start Guide

## iOS Install:

1. Visit the App Store and search for the app “Wi-Viewer HD Pro”
2. Install “Wi-Viewer HD Pro”


## Android Install:

1. Visit the Google Play Market and search for the app “Wi-Viewer HD Pro”
2. Install “Wi-Viewer HD Pro”

**NOTE: Please charge batteries for eight hours before the first use.**

- A. Open the battery door on the SmartMicroScope iGo2 and make note of the SSID.
- B. Turn the SmartMicroScope iGo2 on and leave on for about 10 seconds.
- C. Turn the tablet on and go to SETTINGS and choose WiFi.
- D. Choose the network SSID from the inside battery cover.
- E. Password is “12345678”
- F. After connected, go to the HOME screen and open the application.

\*\*\*\*\* iOS devices ONLY: Touch the ON button on the bottom left of the screen to get the live image.

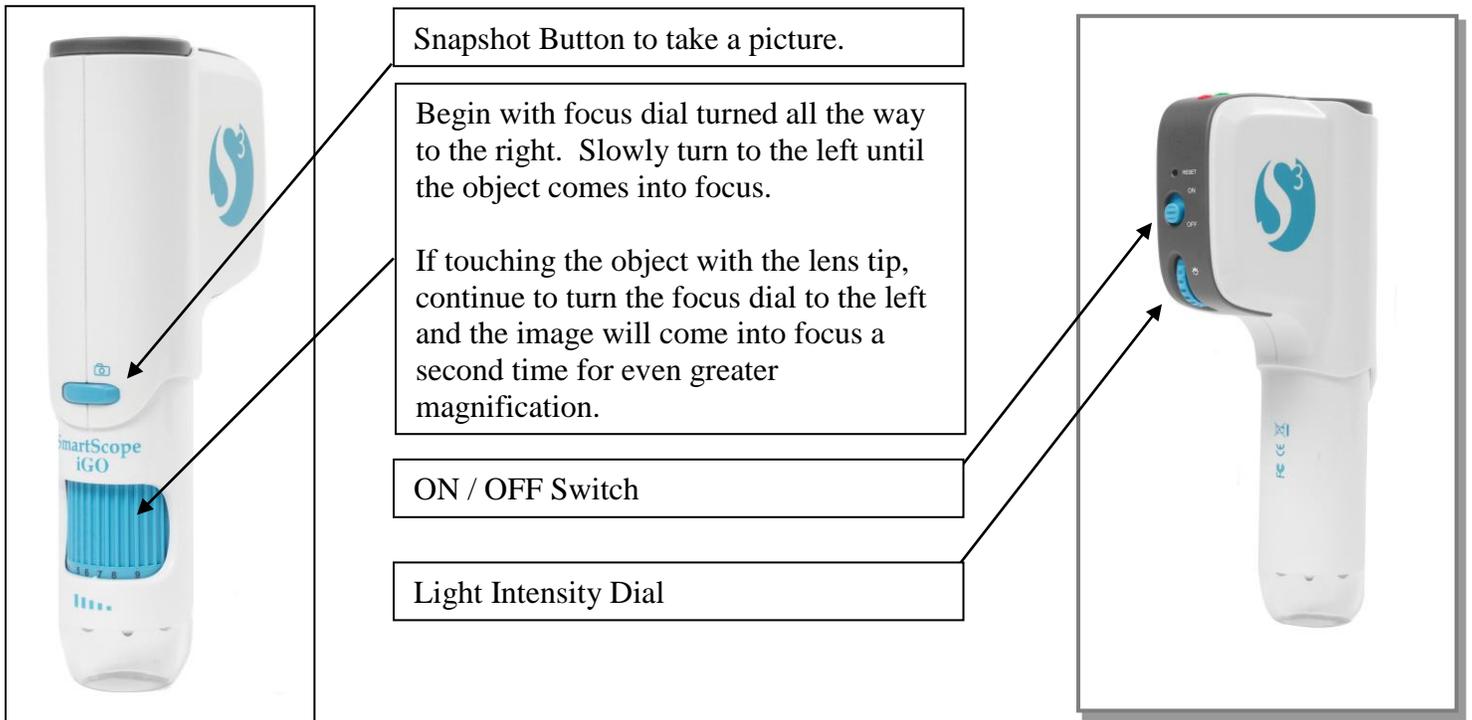

\*\*\*\*\*Technical support, please call 281-312-1297, 1-866-722-2051 or email [support@smartschoolsystems.com](mailto:support@smartschoolsystems.com)